

Turning on the Autoclave



★ Fill Water

If this icon appears please fill water.



⚠ Use mineral free water only.



★ Fill Water

During the filling make sure that the water **does not reach the Red Area**.



Temperature Sensors

Temperature sensors must be inside the bottles.

In case there is only one bottle, one sensor will be inside the bottle and one sensor will be inside the chamber.



Temperature Sensors

Temperature sensors must be inside the bottles.

In case there is more than one bottle, place the two temperature sensors into two separate liquid containers.



For proper sterilization, fill the bottles with approximately the same amount of liquid.

Preparing the Autoclave for Operation


Select the program using the “UP” or “DOWN” arrows.

(Programs can only be selected when the door is open).



Door Closure

Hold the handle and push

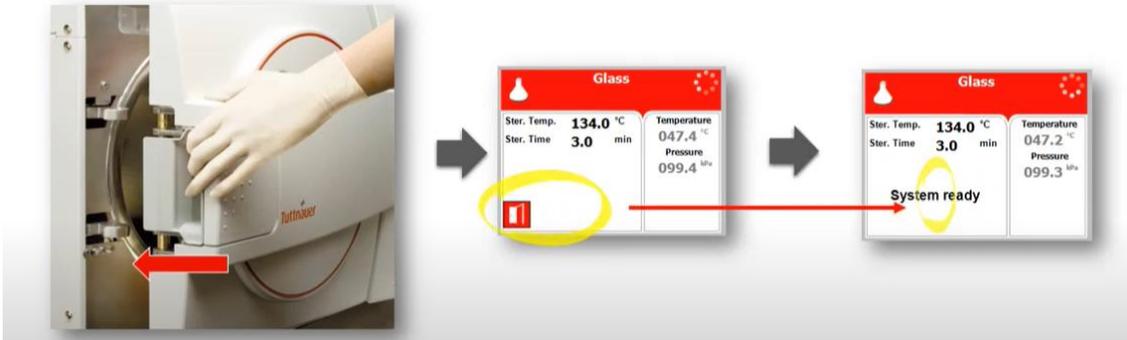


Start Cycle

Press the **START/STOP** key to start the cycle



Cycle

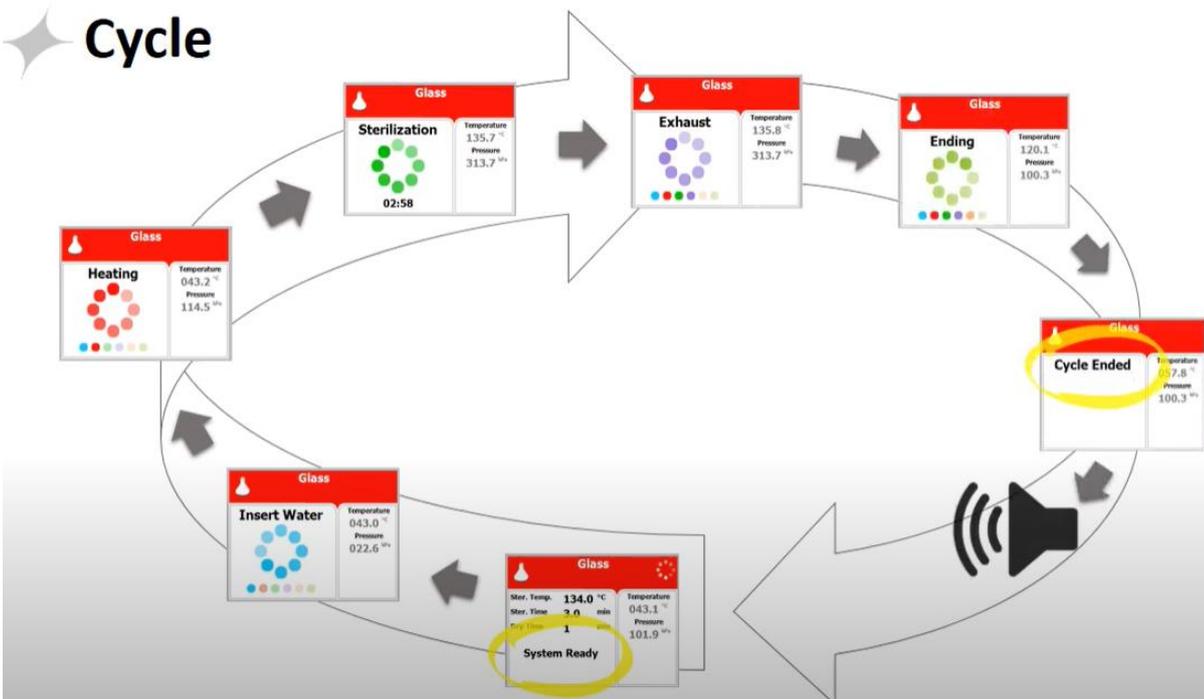

End of Cycle



Press **START/STOP**
key once to confirm

Open the Door



Caution!
The bottles might be hot.
Use heat – resistant gloves